

# Single-assay mRNA vaccine cQA testing based on nanopore sequencing

Steven Verbruggen, Arne Godderis, Anna Matsaroki, Joannes van Cann, Hendrik Van de Voorde, Gerben Menschaert  
OHMX.bio, Noorwegenstraat 4, 9940 Evergem, Belgium, info@ohmx.bio

Strict European guidelines (ICH-Q5 – EMA) regulate QC of biologics as mRNA vaccines. Our extensive service portfolio for mRNA-based therapeutics and vaccines, adeno-associated and other viral vectors, plasmids, microbiomes and/or cell lines, includes analytical solutions based on long-read Nanopore sequencing (Oxford Nanopore Technology, ONT).

We offer a set of assays to assess the combined read-out for several different critical quality attributes (cQAs) of an mRNA drug substance (DS) and drug product (DP).

## mRNA cQAs


### IDENTITY

- Sequence confirmation
- Sequence modifications (m1Ψ and Ψ)

### INTEGRITY

- Intact transcripts
- 5' cap detection
- Poly(A) tail length
- off-target contaminants (as a proxy to dsRNA)

## Sequencing assay overview

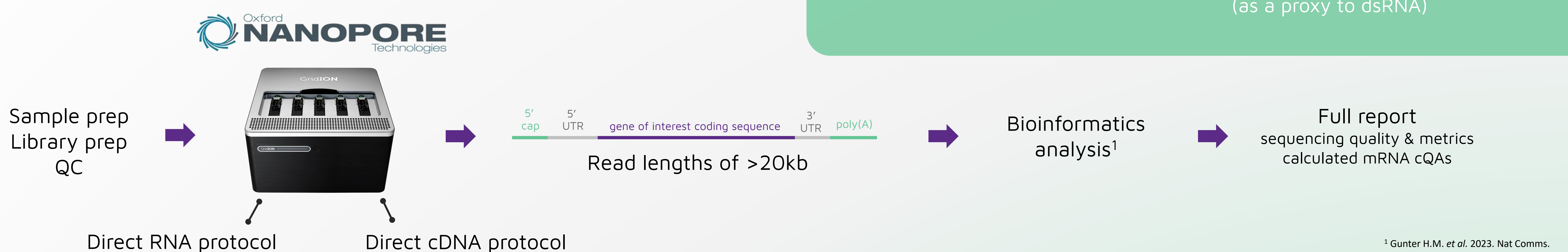


## Full sequence confirmation


- Using Direct cDNA sequencing protocol



### Full mRNA sequence confirmation

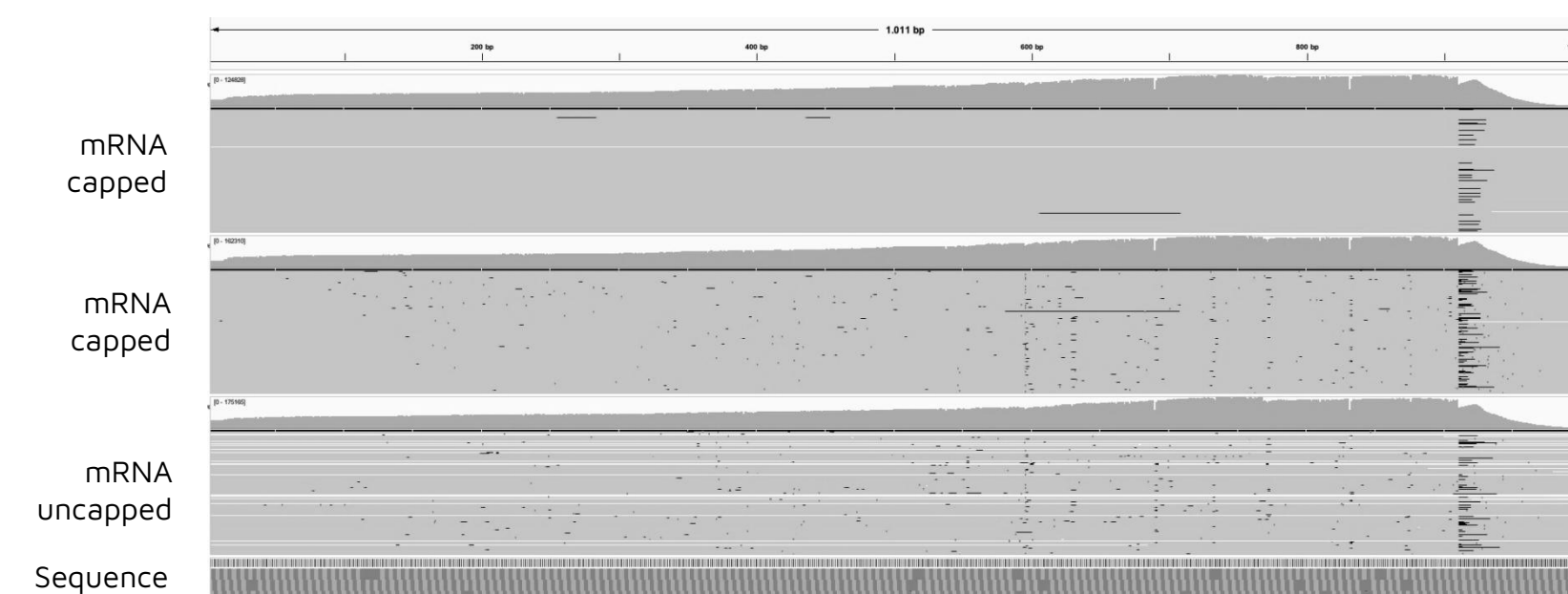


### Sequence length distribution

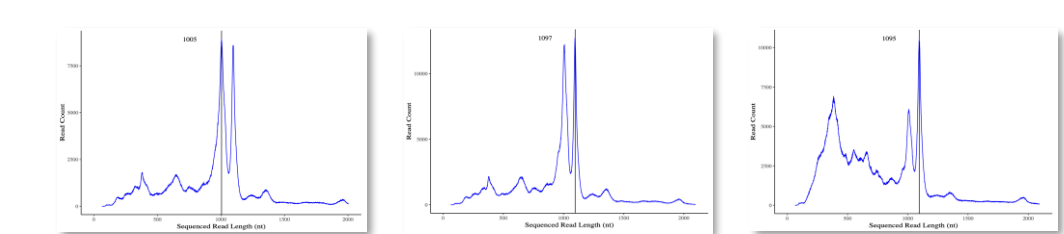


Capping	Reference length	Identified read lengths
Yes	1010	1005 ±1100
Yes	1010	1097 ±1010
No	1010	1095 ±1010

### Full mRNA sequence

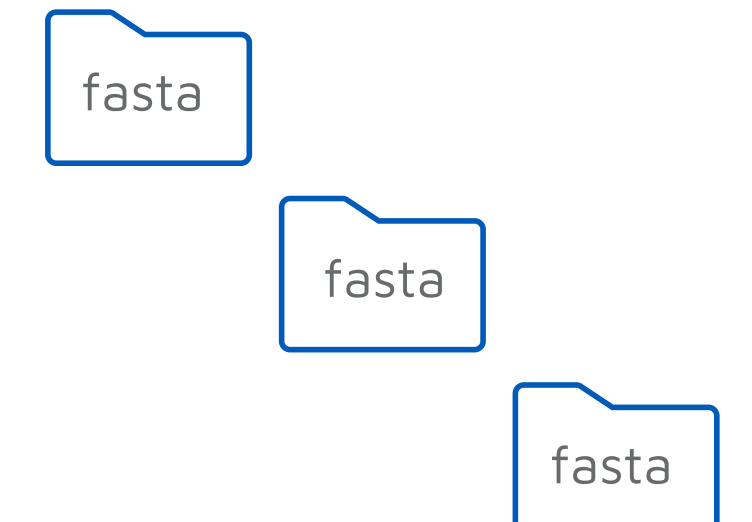


## Sequence modification


- Using Direct RNA sequencing protocol
- Based on deep learning basecalling models specific for m1Ψ and Ψ

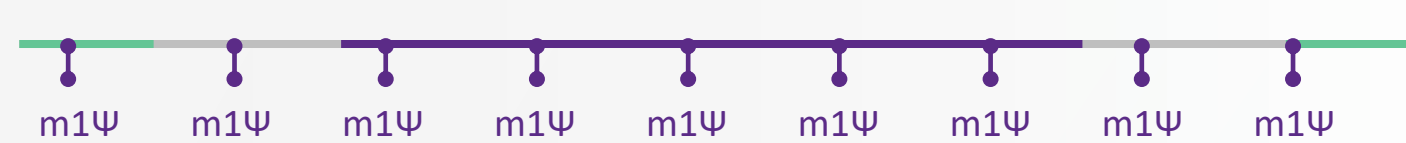

### Full m1-pseudouridine (m1Ψ) modification profile

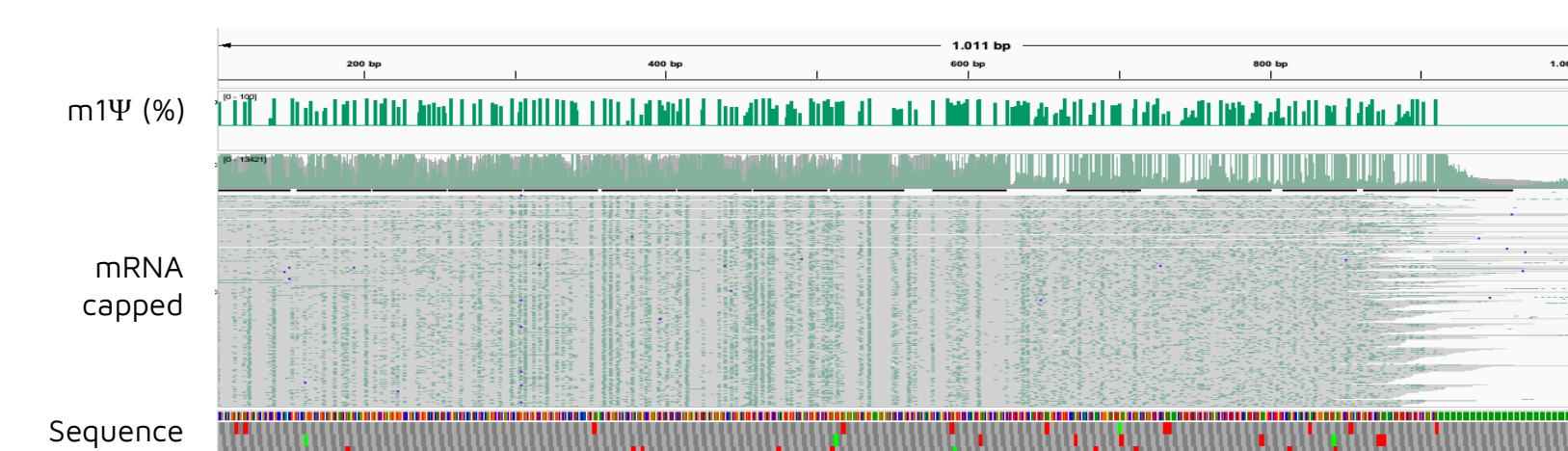


### Detailed zoom in

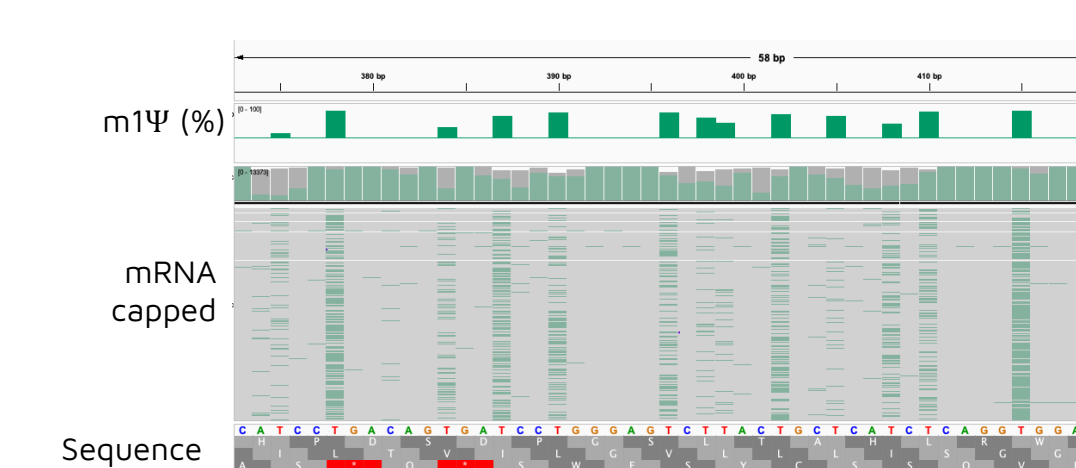


### Modification at base level

Uridine nucleosides in mRNA-vaccines are modified to reduce immunoreactivity and enhance protein production

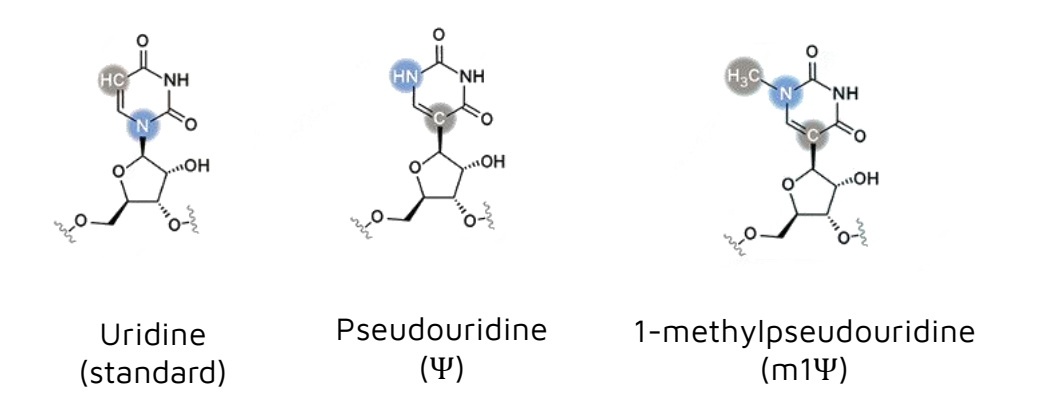


## 3' poly(A) tail / 5' capping


- Using Direct cDNA sequencing protocol
- Based on custom bioinformatics pipelines<sup>2</sup>

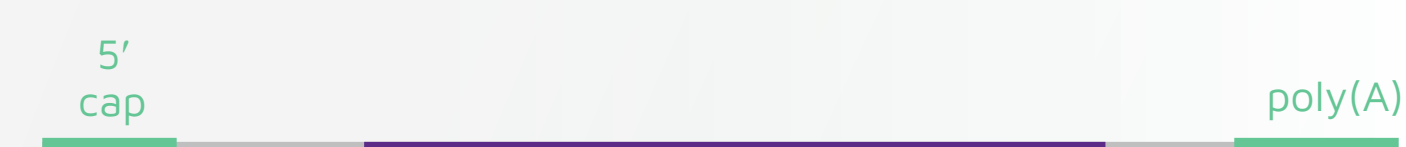

### 5' cap detection

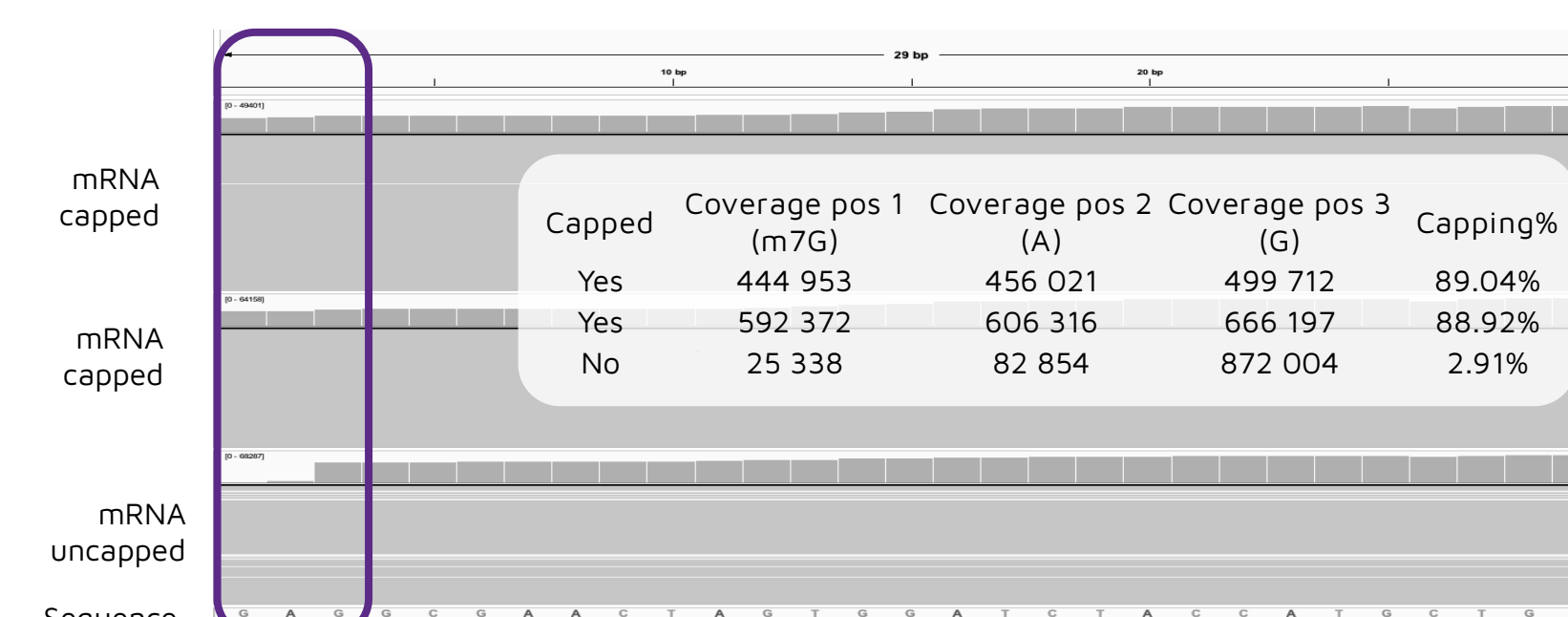


### 3' poly(A) tail

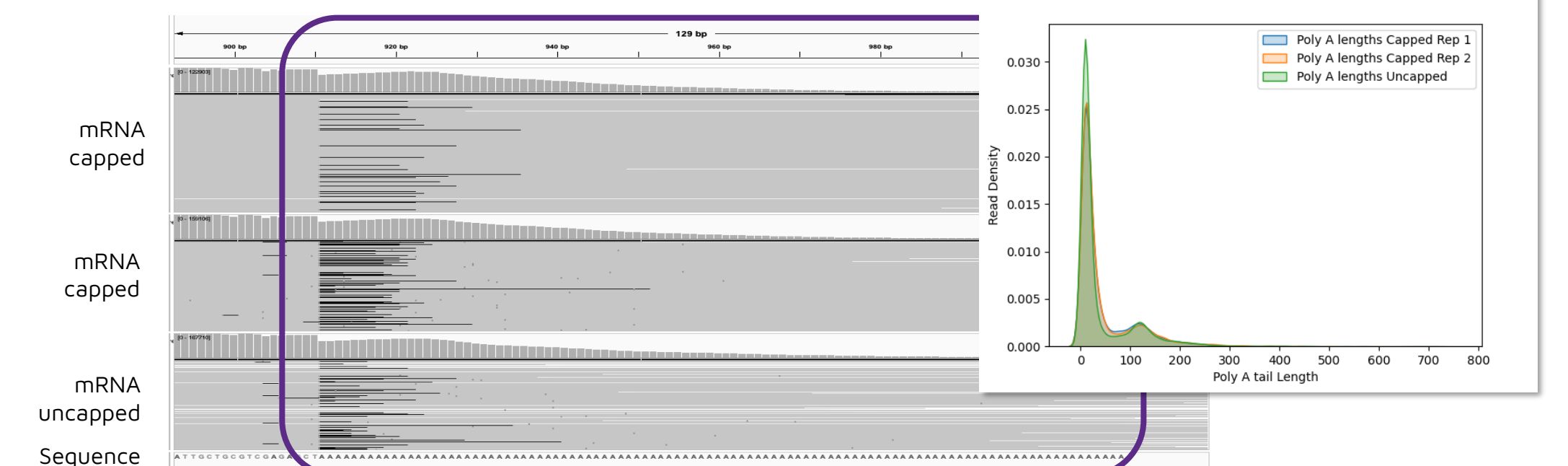


<sup>2</sup> Krause M. et al. 2019. RNA

## Off-target contaminants

- Using Direct RNA sequencing protocol
- Additional poly(A)-tailing
- Based on custom bioinformatics pipelines

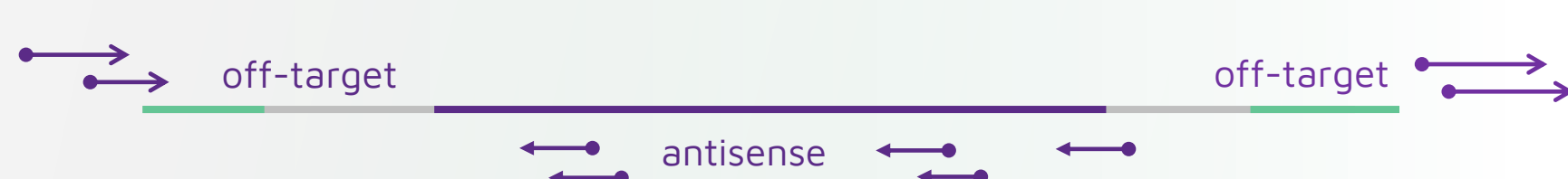

### mRNA species identifications<sup>1</sup>

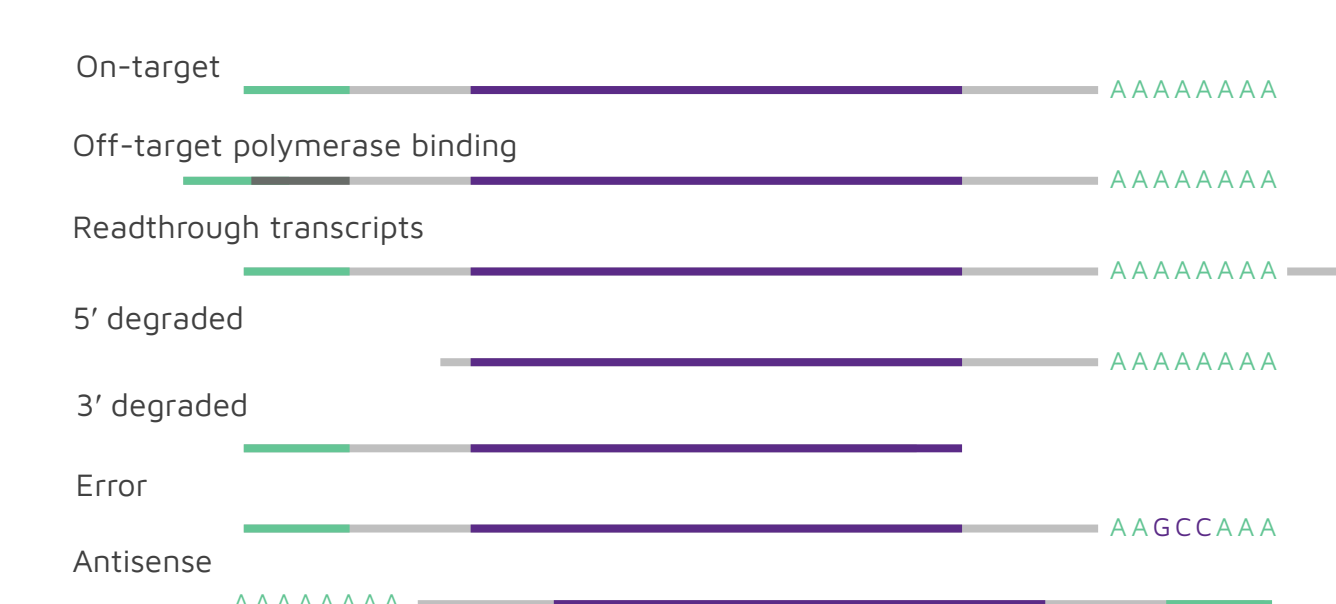

Antisense as proxy to dsRNA %

### On-target and E. coli contaminant mapping

Capped	On-target mapping (%)	E. coli mapping (%)
Yes	99.13%	0.43%
Yes	99.18%	0.27%
No	95.74%	2.94%

<sup>1</sup> Gunter H.M. et al. 2023. Nat Comms

## Towards GMP Oxford Nanopore Q-line GridION system

- Proven, locked-down technology
- Compliance with 21 CFR Part 11 and EU GMP Annex 11 requirements
- Full operating software and consumable version support and supply for three years post-release
- ISO 9001:2015 certified product manufacturing process



- RUO only
- TAT 1-2 weeks
- Non-standardized reporting



- GMP
- TAT <1 week
- Standardized reporting